Equipment Training Sheet - TA.XTPlus Texture Analyser

	Equipment Truming Sheet 171.XTT las Texture Thialysel
Name:	Position:
Department:	Phone:
e-mail:	Director:
Staff Trainee Initials Initials	
	Anyone needing access to MSE equipment needs to contact MSE tech staff: • For initial training • For subsequent usage (will receive equipment card) • No secondary training of the equipment Review of hazards, controls, personal protective equipment:
	 Always wear safety glasses when operating the MTS, also be aware of anyone close by to either move to a safe distance or to wear safety glasses Always use machine shield when testing samples in the MTS Pinch and crush hazard due to moving crosshead, extra caution should be taken Be mindful of the location of the emergency stop button
	 Review of standard operating procedure: http://www.mse.engin.umich.edu/internal/procedures/ta-xtplus-texture-analyser Remember to set the hardware limits on the testing frame to prevent crashing the grips and causing damage to the crosshead and\or load cell Before testing calculate the maximum load conditions of your sample to determine if the machine\load cell is appropriate There is currently a 30kg load cell on the texture analyzer
	 Available supplies to help in the procedure: Rulers and calipers Tweezers Safety glasses Limited amount of test fixtures are available for this tester
Staff Signature/d	ate:

Trainee Signature/date: _____